



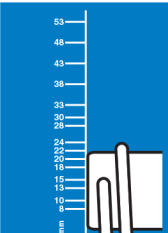

## CAP AND BOTTLE MEASURER

### TO MEASURE CAP:

Place rule across underside of cap with notch flush against inner wall. Read to opposite inner wall of cap.

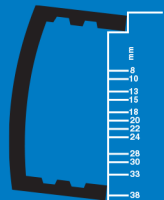

### TO MEASURE BOTTLE NECK FINISH:

Place rule across finish with measurer tightly against thread. Read to outer edge of thread on opposite side of finish.



A Division of Berlin Packaging

Corporate One West  
1195 Washington Pike  
Bridgeville, PA 15017-2854

**1.800.922.7558**

412.257.3100

F 412.257.3001

qorpak.com

An ISO 9001:2008  
Certified Company

SEE OTHER SIDE



## COMMON SIZE INFORMATION

Screw thread closure sizes are expressed in two numbers.

The first number refers to the closure diameter (in mm) and the second number represents the GPI finish (the top of the container).

For example, a neck size of **28-430** has a diameter of **28 mm** and fits a container with a **430 GPI thread finish**.



## COMMON CAP FINISHES

